

PB.5047

scale magnifier



euromex
microscopes holland

user
manual


Thank you for purchasing the PB.5047 scale magnifier. It is designed to have a wide angle, high imaging definition and high measurement accuracy

Instructions

- 1.** By rotating the focus adjustment ring (A), the scale and the object to be measured can be put in to focus
- 2.** If needed the LED illumination can be turned on and off by switching the white button (C)

To change the batteries:

- A.** Turn the focus adjustment ring (A) to the right, all the way until the two parts of the magnifier are separated (A and B)
- B.** Change the batteries (4pcs of AG9 1.5v)

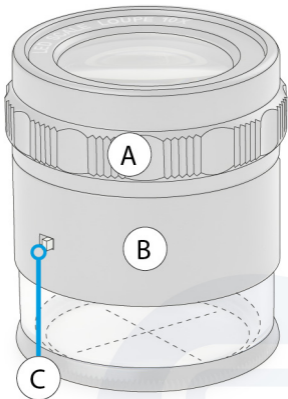

v.425501

*All information may be
changed without prior notice*



