



motorized models



supplementary
user manual



v.424361

Additional instruction for motorized models of Delphi-X

Control buttons

- A. Objective selection buttons
- B. Function button
- C. Side buttons

Operation

There are two switching modes:

1. Free switching mode

In the free switching mode, all switching buttons (A) are available for use. The buttons are lit green. Pressing the predefined buttons on the front of the Delphi-X Observer microscope will select the corresponding objective. By default, from left to right, the order is: 4x/10x/20x/40x/100x. When using the side buttons (C) all objectives can be chosen clockwise and counterclockwise

2. Fast switching mode (switching between two selected objectives)

In the fast switching mode only the buttons of the selected objectives are available for use

- Green (active objective)
- Yellow (the other selected objective)


When using the side buttons (C) these two objectives can be chosen

Selecting the two objectives for switching mode:

- Press and hold the green function button (B). Now all indicator lights of switching buttons will be turned off
- Then press the two buttons of the preferred objectives in the preferred order. Those two indicator lights will turn on (one green, one yellow)
- Finally, release the "function button"

The side buttons (C) can be used to, either switch between all objectives, or the two selected objectives, depending on the selected mode. The function of the left side and right side buttons are the same



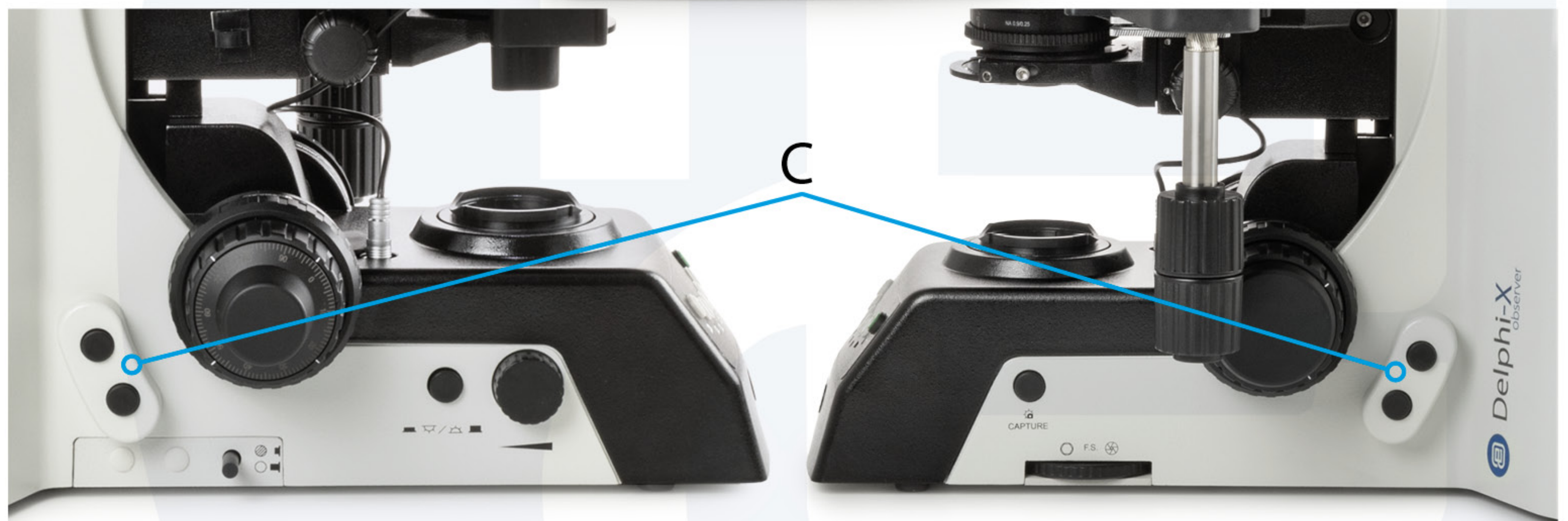

Euromex Microscopen bv • Typograaf 8 • 6921 VB Duiven • The Netherlands
 T +31 (0) 26 323 22 11 • info@euromex.com • www.euromex.com



All information may be changed without prior notice

