DC.5000-WiFi-2



Contents

The DC.5000-WiFi-2	2
Connecting to the camera's WiFi	2
Connect to WiFi with Windows 10	2
Connect to WiFi with Windows 7/8	3
Connect to WiFi with MAC os	4

The DC.5000-WiFi-2

This camera can run in either:

- USB-2 mode
- WiFi mode

The mode can be selected by pushing the **MODE** button

<u>Please note:</u> the camera can not run in both ways simultaneously. The USB indicator LED will light up if the camera is set in USB-2 mode and the WiFi indicator LED will light up when set in WiFi mode

 In USB-2 mode: connect the USB cable to the PWR/USB connection on the camera and connect the other side to your computer.
 The camera now functions like all Euromex USB cameras and uses ImageFocus alpha.
 See dedicated ImageFocus alpha manual for operation in this mode



In WiFi mode: connect the USB cable to the PWR/USB connection on the camera and connect the other side
to the power adapter. The camera now functions like a WiFi camera and uses ImageFocus alpha PC software or
the ImageFocus alpha app's

Connecting to the camera's WiFi

The first step that needs to be taken on any device is connecting to the WiFi camera's own network. This means you will need to disconnect from your standard WiFi network. The instructions for this are shown below

After connection to the camera's WiFi network please refer to the dedicated ImageFocus alpha manual for use with PC, and decicated manuals for the app's. The IOS app can be downloaded from the app store, for Android a download is available on the www.euromex.com download section

Connect to WiFi with Windows 10

 On your desktop, click the network icon in bottom right-hand corner of your screen. This will bring up a list of available connections



- 2. Find the name (SSID) of the CMEX WiFi network and click on it
- 3. Ensure connect automatically is checked and then click connect
- 4. Enter your WiFi password. Standard password is: 12345678
- 5. Click OK to finish. Your computer should remember your password for you from now on

Connect to WiFi with Windows 7/8


1. Open the desktop



Click the network icon in the bottom right-hand corner of your screen. This will bring up a list of available network connections



3. Find the name (SSID) of your CMEX WiFi network and click on it

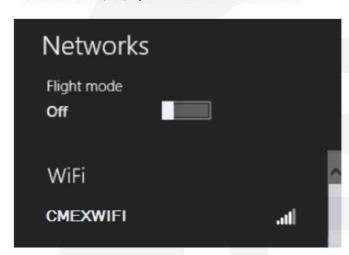

4. Ensure connect automatically is checked and then click the connect button.


Enter your WiFi password, and then click OK to finish. Your computer should remember your password for you from now on. Standard password is: 12345678



Connect to WiFi with MAC os

Connect using the AirPort®/WiFi menu

1. On the desktop, click the AirPort/WiFi icon - - -, then select the CMEX WiFi name (SSID) you want to connect to

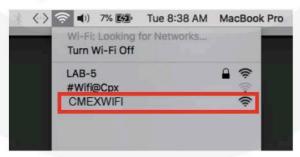

2. When required, enter the CMEX WiFi password then click join. Standard password is: 12345678



Euromex Microscopen bv • Papenkamp 20 • 6836 BD Arnhem • The Netherlands T +31 (0) 26 323 22 11 • info@euromex.com • www.euromex.com







