

Euromex cameras



1. Software installation

In order to use the Euromex cameras, you need to install the application software, available and downloadable (no charge) on the Euromex website (downloads).

Differences in function and appearance between Windows and Mac apply

1.1 Windows

- Select and download the ImageFocus software for Windows


Important:

When you download the software from the Euromex website, please first unpack the file to a location on your harddrive first(not on the desktop!), then run the installation file

- Right mouse click and run the file with "Run as administrator". If running as administrator is not possible on your computer, please run the file by left mouse clicking the file

1.2 Mac

- Select and download ImageFocus software for Mac
- Double-click on the zip file to unpack. double-click on the dmg-file
- Drag IFA icon to your applications folder

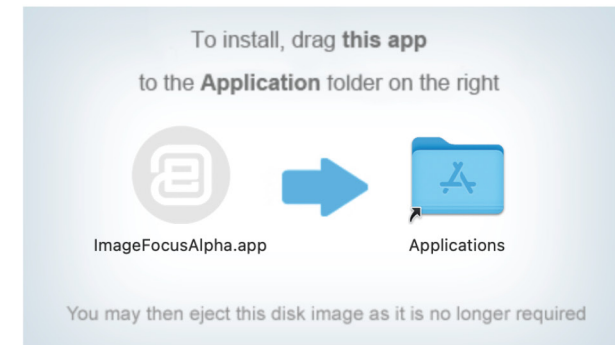


1.3 How to start the camera


ImageFocus will detect all of the cameras that your computer has installed and will list all the camera names. Click on the camera name to run the camera

1.4 Full manual (English only) available in the software:

- ImageFocus Plus

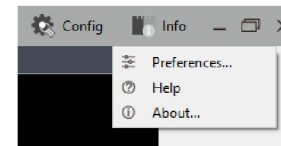

- ImageFocus 4

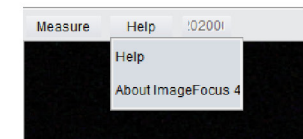



EN: Standard c-mount adapter for tube 23.2 mm
 DE: 23,2 mm Tubus Standard C-mount Adapter
 FR: Adaptateur de montage en C standard pour tube de 23,2 mm
 ES: Adaptador estándar de montura c para tubo de 23,2 mm
 IT: Adattatore standard per attacco a C per tubo da 23,2 mm
 NL: Standaard c-mount adapter voor tubus 23,2 mm



EN: Model specific c-mount adapter
 DE: Modellspezifischer C-mount Adapter
 FR: Adaptateur de montage en C spécifique au modèle
 ES: Adaptador de montaje en c específico del modelo
 IT: Adattatore per montaggio a C specifico per il modello
 NL: Modelspecifieke c-mount adapter



EN: Brand specific c-mount adapter
 DE: Markenspezifischer C-mount Adapter
 FR: Adaptateur de montage en C spécifique à la marque
 ES: Adaptador de montaje en c específico de la marca
 IT: Adattatore per montaggio a C specifico per la marca
 NL: Merkspecifieke c-mount adapter

23.2 mm tubes

30 / 30.5 mm tubes

